

Anti-Laminin-511, Human, Mouse-Mono (Clone 12D)

Catalog NO. FDV-0026

Research use only, not for human or animal therapeutic or diagnostic use.

Product Background

Laminins, which consist of three subunits called α , β and γ chains, are major cell-adhesive components of extracellular matrix, especially basement membranes (BMs). The laminin family is constituted of over 15 isoforms, and each member is expressed in a tissue-specific manner and plays a differential role in each tissue. Laminin-511 (Lm511) (formerly laminin-10), which consists of $\alpha 5$, $\beta 1$ and $\gamma 1$ chains, appears at an early stage of development and is most widely expressed in adult tissues, mainly basement membranes of various epithelial tissues and vasculatures. Recently, laminin-511 is often used as a culture substrate for ES and iPS cells.

A mouse monoclonal antibody (clone 12D) for laminin-511, was established using human placental laminin as an antigen. Clone 12D specifically interacts with the trimeric structure of laminin-511 isoform but does not react with any of the $\alpha 5$, $\beta 1$ and $\gamma 1$ chains in the reducing conditions. Furthermore, it has been validated that clone 12D does not react with any of laminins-111, -211, -332, -3B32, -411 and -521. Unlike conventional anti- $\alpha 5$ antibodies, clone 12D distinguishes laminin-511 from laminin-521. This is an important tool to evaluate specific distribution and physiological functions of laminin-511. In human cancers, clone 12D detects small vascular vessels with high sensitivity and hence useful for evaluating tumor angiogenesis.

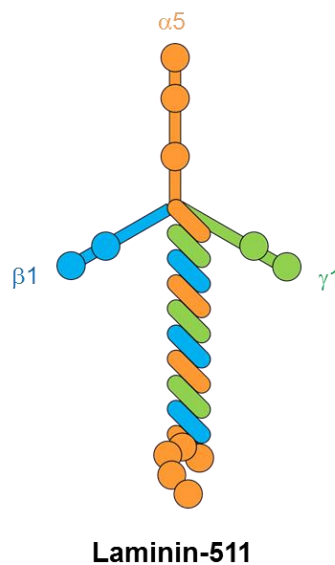


Figure 1. Protein structure of laminin-511

Description

Catalog Number : FDV-0026

Format : Mouse ascites

Volume : 100 μ L

Formulation : Ascites without any additives

Host Species and Clonality : Mouse monoclonal

Isotype and Subclass : IgG2b

Lot Number : see vial label

Specificity : Human, other species not tested yet

Storage : For short-term storage, -20°C. For long-term storage, -80°C storage is preferable.

Avoid repeated freeze-thaw cycles and avoid storage at 4°C.

Application

- Western blotting under non-reducing conditions
 - Immunohistochemistry under frozen section
 - Immunoprecipitation and immunoaffinity purification
 - ELISA
-

Recommended usage

- Western blotting 1/1,000
 - Immunohistochemistry Optimal dilutions should be determined by the researcher.
 - Immunoprecipitation Optimal conditions should be determined by the researcher.
 - ELISA Optimal conditions should be determined by the researcher.
-

Application examples


Western blot

Sample : Recombinant human intact laminin-511 (Lm511)
consisting of α 5, β 1 and γ 1 chains

Gel conc. : 4.5% under non-reducing conditions


1st Antibody : clone 12D, 1/1,000 dilution

2nd Antibody : anti-mouse IgG (H+L)-HRP conjugate

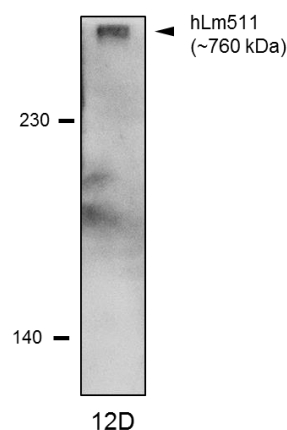

Specificity of 12D for laminin isoforms

Seven recombinant human laminins were coated on a multi-well plate and detected by clone 12D. High specificity of clone 12D to Lm511 is observed.

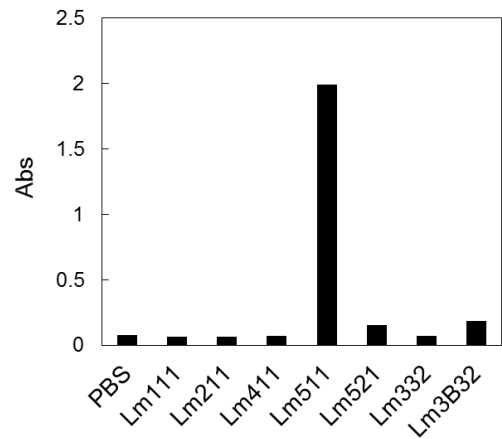


Immunohistochemistry : Human normal mammary gland

Sample : Human mammary gland (frozen slice)

Black arrows : Vascular basement membrane

White arrowheads : Mammary gland basement membrane

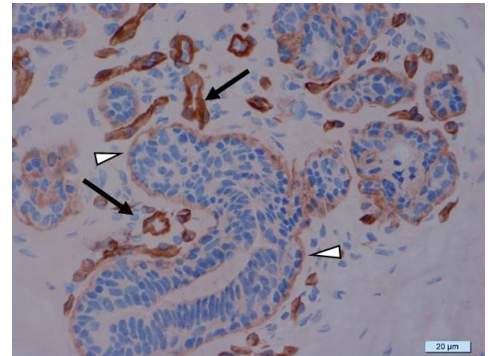


Immunohistochemistry : Breast cancer tissue

Sample : Non-invasive carcinoma tissue (frozen slice)

A: Vascular basement membrane


B: Mammary gland basement membrane

C: Basement membrane-like structures surrounding tumor cells

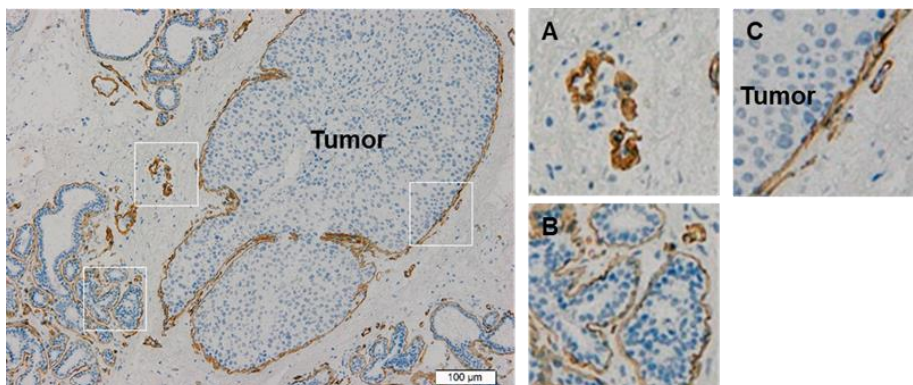

Immunohistochemistry : Lung cancers

Sample : Lung squamous cell carcinoma (left) and lung adenocarcinoma (right)

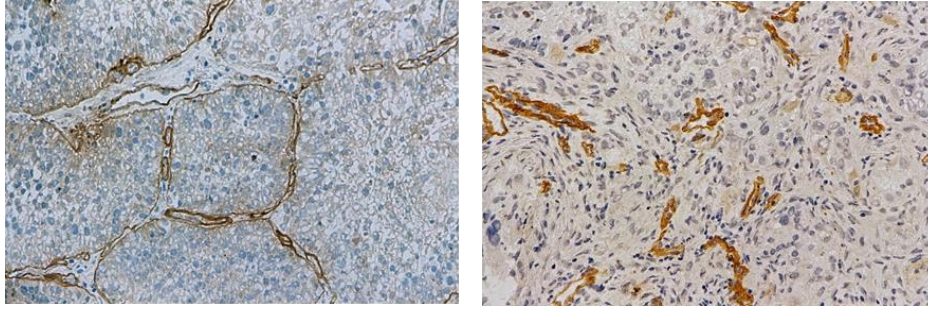

*All data are provided from Dr. Kaoru Miyazaki

Reference

1. Komiya *et al.*, *Cancer Med.*, **3**, 537-549 (2014) Angiomodulin (AGM/IGFBP-rP1), a marker of cancer vasculature, is upregulated by vascular endothelial cell growth factor and increases vascular permeability as a ligand of integrin $\alpha v \beta 3$.

Related products

Catalog No.	Product name	Target	Application
FDV-0023	Anti-Laminin $\alpha 3B$, Human, Mouse-Mono (F7)	Laminin $\alpha 3B$	IHC, WB, IP, ELISA
FDV-0024	Anti-Laminin $\alpha 3A$, Human, Mouse-Mono (BG5)	Laminin $\alpha 3A$	IHC, WB, IP, ELISA
FDV-0025	Anti-Laminin $\gamma 2$ N-terminal fragment, Human, Mouse-Mono (P2H)	Laminin $\gamma 2$ N-terminal fragment	IHC, WB, ELISA
FDV-0026	Anti-Laminin 511, Human, Mouse-Mono (12D)	Trimeric Lm511 structure	IHC, WB, IP, ELISA

Disclaimer/免責事項

This product has been commercialized by Funakoshi Co., Ltd. based on the results of academic research, and the advertisement text, figures and manuals (hereinafter "Product information") have been prepared based on published research reports on December, 2018. The academic interpretation at the time of creation of the Product Information may change in accordance with future developments in the relevant research field and expansion of various scientific findings, and the latest version and certainty of the Product Information are not guaranteed. The specifications of this product and the Product Information are subject to change without notice. Please contact us for the latest information.

本商品は学術研究成果を基にフナコシ株式会社が商品化したもので、2018年12月時点における公開研究報告を基に広告文章およびマニュアル(以下、商品資料)を作成しています。今後の当該研究分野の発展および各種学術知見の拡大にともない、商品資料作成時の学術的解釈が変更になる可能性があり、最新性・確実性を保証するものではありません。また、本商品の仕様および商品資料を予告なく変更する場合がございます。最新の情報に関しましては、弊社までご確認いただけますようお願い申し上げます。

 E-mail Newsletter Sign Up

Japanese 

English 

 **funakoshi**
FRONTIERS IN LIFE SCIENCE

URL: <http://funakoshi.co.jp>
9-7-1 Hongo 2-Chome, Bunkyo-ku, Tokyo 113-0033