Product Manual

CytoSelect™ 24-Well Cell Invasion Assay (Collagen I, Fluorometric Format)

Catalog Number

CBA-111-COL 12 assays

FOR RESEARCH USE ONLY Not for use in diagnostic procedures




Introduction

The ability of malignant tumor cells to invade normal surrounding tissue contributes in large part to the significant morbidity and mortality of cancers. Invasiveness requires several distinct cellular functions including adhesion, motility, detachment, and extracellular matrix proteolysis. Metastatic cells produce many proteolytic enzymes (e.g. lysosomal hydrolysates, collagenases, plasminogen activators) while the expression of certain cell surface protease receptors is also increased.

Cell Biolabs CytoSelectTM Collagen Cell Invasion Assay Kit utilizes Bovine Type I Collagen-coated inserts to assay the invasive properties of tumor cells. It contains sufficient reagents for the evaluation of 12 samples.

Assay Principle

The CytoSelectTM Collagen Cell Invasion Assay Kit contains polycarbonate membrane inserts (8 µm pore size) in a 24-well plate. The upper surface of the insert membrane is coated with a uniform layer of dried Bovine Type I Collagen matrix. This collagen matrix layer serves as a barrier to discriminate invasive cells from non-invasive cells. Invasive cells are able to degrade the collagen matrix layer, and ultimately pass through the pores of the polycarbonate membrane. Finally, these invaded cells are then dissociated from the membrane and subsequently detected by the patented CyQuant® GR Dye (Invitrogen).

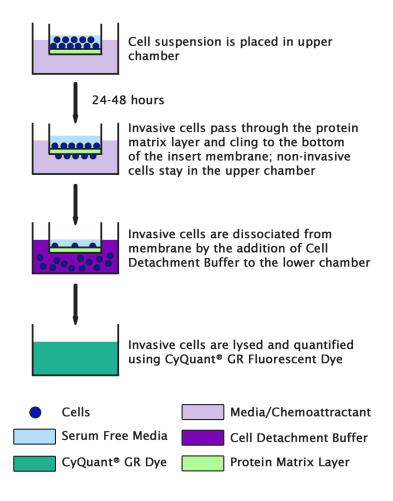



Related Products

- 1. CBA-101-C: CytoSelect™ 24- Well Cell Migration and Invasion Assay (8µm, Fluorometric)
- 2. CBA-106-C: CytoSelectTM 96-Well Cell Migration and Invasion Assay (8μm, Fluorometric)
- 3. CBA-111: CytoSelectTM 24-Well Cell Invasion Assay (Basement Membrane, Fluorometric)
- 4. CBA-112: CytoSelectTM 96-Well Cell Invasion Assay (Basement Membrane, Fluorometric)
- 5. CBA-112-COL: CytoSelectTM 96-Well Cell Invasion Assay (Collagen I, Fluorometric)

Kit Components

- 1. <u>Collagen Invasion Chamber Plate</u> (Part No. 111001-COL): One 24-well plate containing 12 collagen-coated cell culture inserts
- 2. <u>Cell Detachment Solution</u> (Part No. 10101): One 5 mL bottle
- 3. 4X Lysis Buffer (Part No. 10102): One 5 mL bottle
- 4. CyQuant® GR Dye (Part No. 10103): One 25 μL tube
- 5. Forceps (Part No. 11005): One each

Materials Not Supplied

- 1. Invasive cell lines
- 2. Cell culture medium
- 3. Serum free medium, such as DMEM containing 0.5% BSA, 2 mM CaCl₂ and 2 mM MgCl₂
- 4. Cell culture incubator (37°C, 5% CO₂ atmosphere)
- 5. 96-well plate suitable for a fluorescence plate reader
- 6. Fluorescence plate reader

Storage

Store all components at 4°C.

Assay Protocol

- 1. Under sterile conditions, allow the collagen invasion chamber plate to warm up at room temperature for 10 minutes.
- 2. Rehydrate the collagen layer of the cell culture inserts by adding 300 μ L of warm, serum-free media to the inner compartment. Incubate at room temperature for 30 minutes.
- 3. Prepare a cell suspension containing $0.5-1.0 \times 10^6$ cells/ml in serum free media. Agents that inhibit or stimulate cell invasion can be added directly to the cell suspension.
 - *Note: Overnight starvation may be performed prior to running the assay*
- 4. Carefully remove 250 μ L of rehydration medium (step 2) from the inserts without disturbing the collagen layer (leaving 50 μ L inside).



- 5. Add 250 µL of the cell suspension solution to the inside of each insert.
- 6. Add 500 μL of media containing 10% fetal bovine serum or desired chemoattractant(s) to the lower well of the invasion plate.
- 7. Incubate for 12-24 hours in a cell culture incubator.
- 8. Carefully aspirate the media from the inside of the insert. Transfer the insert to a clean well containing 225 μ L of Cell Detachment Solution. Incubate 30 minutes at 37 °C.
- 9. Completely dislodge the invaded cells from the underside of the membrane by gently tilting the insert several times in the detachment solution. Remove and discard the insert.
- 10. Prepare sufficient 4X Lysis Buffer/CyQuant® GR dye solution for all samples by diluting the dye 1:75 in 4X Lysis Buffer (for example, add 5 μ L to 370 μ L of 4X Lysis Buffer).
- 11. Add 75 µL of 4X Lysis Buffer/CyQuant® GR dye solution to each well containing cells and 225 µL of Cell Detachment Solution. Incubate 20 minutes at room temperature.
- 12. Transfer 200 μ L of the mixture to a 96-well plate suitable for fluorescence measurement. Read fluorescence with a fluorescence plate reader at 480 nm/520 nm.

Example of Results

The following figures demonstrate typical with the CytoSelectTM Collagen Cell Invasion Assay Kit. One should use the data below for reference only. This data should not be used to interpret actual results.

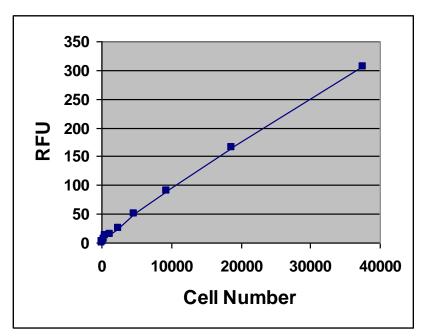


Figure 1. Quantitation of Human HT-1080. HT-1080 cell suspension was titrated 1:2 in Cell Detachment Buffer. 150 μ L of each dilution was transferred to a 96-well, fluorometer plate, followed by addition of 50 μ L 4X Lysis Buffer/CyQuant® GR dye solution (1:75).



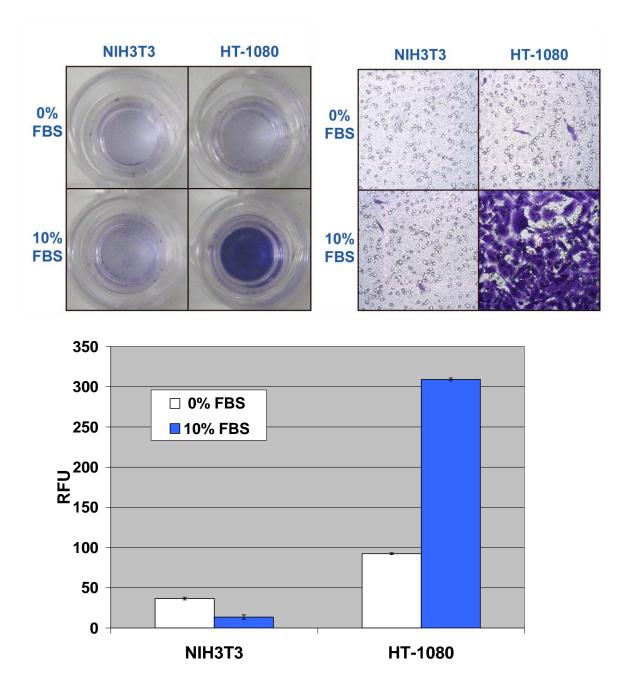


Figure 2. Human Fibrosarcoma HT-1080 Collagen Cell Invasion. HT-1080 and NIH3T3 (negative control) were seeded at 200,000 cells/well and allowed to invade toward FBS for 24 hrs. Invasive cells on the bottom of the invasion membrane were stained (top pictures) and quantified by CyQuant® GR dye as described in the Assay Protocol (bottom figure).

References

- 1. Erkell, L. J., Schirrmacher, V. (1988) Cancer Res 48, 6933-6937.
- 2. Montgomery, A. M. P., De Clerck, Y. A., Langley, K. E., Reisfeld, R. A., Mueller, B. M. (1993) *Cancer Res* **53**,693-700.
- 3. Monsky, W. L., Lin, C. Y., Aoyama, A., Kelly, T., Akiyama, S. K., Mueller, S. C., Chen, W. T. (1994) *Cancer Res* **54**,5702-5710.



Recent Product Citations

- 1. Zhen, Q. et al. (2020). EPAS1 promotes peritoneal carcinomatosis of non-small-cell lung cancer by enhancing mesothelial-mesenchymal transition. *Strahlenther Onkol*. doi: 10.1007/s00066-020-01665-1.
- 2. Negro, G. et al. (2020). Molecular heterogeneity in breast carcinoma cells with increased invasive capacities. *Radiol Oncol.* **54**(1):103-118. doi: 10.2478/raon-2020-0007.

License Information

This product is provided under an intellectual property license from Life Technologies Corporation. The purchase of this product conveys to the buyer the non-transferable right to use the purchased product and components of the product only in research conducted by the buyer (whether the buyer is an academic or for-profit entity). The sale of this product is expressly conditioned on the buyer not using the product or its components, or any materials made using the product or its components, in any activity to generate revenue, which may include, but is not limited to use of the product or its components: (i) in manufacturing; (ii) to provide a service, information, or data in return for payment; (iii) for therapeutic, diagnostic or prophylactic purposes; or (iv) for resale, regardless of whether they are resold for use in research. For information on purchasing a license to this product for purposes other than as described above, contact Life Technologies Corporation, 5791 Van Allen Way, Carlsbad CA 92008 USA or outlicensing@lifetech.com.

Warranty

These products are warranted to perform as described in their labeling and in Cell Biolabs literature when used in accordance with their instructions. THERE ARE NO WARRANTIES THAT EXTEND BEYOND THIS EXPRESSED WARRANTY AND CELL BIOLABS DISCLAIMS ANY IMPLIED WARRANTY OF MERCHANTABILITY OR WARRANTY OF FITNESS FOR PARTICULAR PURPOSE. CELL BIOLABS' sole obligation and purchaser's exclusive remedy for breach of this warranty shall be, at the option of CELL BIOLABS, to repair or replace the products. In no event shall CELL BIOLABS be liable for any proximate, incidental or consequential damages in connection with the products.

Contact Information

Cell Biolabs, Inc. 5628 Copley Drive San Diego, CA 92111

Worldwide: +1 858-271-6500 USA Toll-Free: 1-888-CBL-0505 E-mail: tech@cellbiolabs.com

www.cellbiolabs.com

©2006-2024: Cell Biolabs, Inc. - All rights reserved. No part of these works may be reproduced in any form without permissions in writing.

