CarbPROFILE gel



Description

CarbPROFILE gels are monosaccharides-Sepharose affinity matrices used for purification of specific carbohydrate-binding proteins¹. The carbohydrates are attached through their non reducing hydroxyl group after pre-activation of sepharose matrix by divinylsulfone (DVS) (*see scheme 1 below*). The binding of lectins and carbohydrates binding proteins to carbohydrate affinity gel is non-covalent and reversible with high capacity (over 1 mg/mL). Lectins and carbohydrates binding proteins can be recovered by competitive elution (*i.e.* 0.2 to 0.5 M of monosaccharide) or by modulations of pH and/or ionic strength in high yield and purity. Each new batch of affinity matrix produced are fully characterized by measurement of its capability to bind lectins and recover the same amount of proteins from batch to batch which assure a total reliability.



Scheme 1. CarbPROFILE gel matrix

Specifications of CarbPROFILE gel

See below, Table 1, the specification of CarbPROFILE gel matrix.

Reference	Name	Specificity	Capacity (mg of protein/mL of gel)	Unit sizeª <i>(mL)</i>
CGF	Fucose-CarbPROFILE gel	Fucose binding protein	> 15 (based on UEA-I lectin)	5, 10 or 25
CGGa	Galactose-CarbPROFILE gel	Galactose binding protein	> 15 (based on AIA lectin)	5, 10 or 25
CGGN	N-Acetylglucosamine- CarbPROFILE gel	N-Acetylglucosamine binding protein	> 15 (based on WGA lectin)	5, 10 or 25
CGM	Mannose-CarbPROFILE gel	Mannose binding protein	> 30 (based on ConA lectin)	5, 10 or 25
CGR	Rhamnose-CarbPROFILE gel	Rhamnose binding protein	> 15 (based on CorM lectin)	5, 10 or 25

Table 1. Specifications of CarbPROFILE gel. a. available in a pre-packed column or in suspension.

Applications

CarbPROFILE gel is used for purification of specific carbohydrate-binding proteins from naturals or recombinant sources:

- lectins,
- lectins-like,
- GAG-binding proteins.

Distributed by: CliniSciences Group

References

1. Andon, N. L., Eckert, D., Yates III, J. R., Haynes, P. A. Proteomics, 2003, 3, 1270-1278.

CliniSciences Group

Austria

Company: CliniSciences GmbH Address: Sternwartestrasse 76, A-1180 Wien - Austria Telephone: +43 720 115 580 Fax: +43 720 115 577 Email: oesterreich@clinisciences.com Web: https://www.clinisciences.com

Finland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: suomi@clinisciences.com Web: https://www.clinisciences.com

Iceland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: island@clinisciences.com Web: https://www.clinisciences.com

Netherlands

Company: CliniSciences B.V. Address: Kraijenhoffstraat 137A, 1018RG Amsterdam, - Netherlands Telephone: +31 85 2082 351 Fax: +31 85 2082 353 Email: nederland@clinisciences.com Web: https://www.clinisciences.com

Portugal

Company: Quimigen Unipessoal LDA Address: Rua Almada Negreiros, Lote 5, Loja 14, 2615-275 Alverca Do Ribatejo - Portugal Telephone: +351 30 8808 050 Fax: +351 30 8808 052 Email: info@quimigen.com Web: https://www.quimigen.pt

Switzerland

Company: CliniSciences AG Address: Fracht Ost Flughafen Kloten CH-8058 Zürich - Switzerland Telephone: +41 (044) 805 76 81 Fax: +41 (044) 805 76 75 Email: switzerland@clinisciences.com Web: https://www.clinisciences.com

Belgium

Company: CliniSciences S.R.L Address: Avenue Stalingrad 52, 1000 Brussels - Belgium Telephone: +32 2 31 50 800 Fax: +32 2 31 50 801 Email: belgium@clinisciences.com Web: https://www.clinisciences.com

France

Company: CliniSciences S.A.S Address: 74 Rue des Suisses, 92000 Nanterre- France Telephone: +33 9 77 40 09 09 Fax: +33 9 77 40 10 11 Email: info@clinisciences.com Web: https://www.clinisciences.com

Ireland

Company: CliniSciences Limited Address: Ground Floor, 71 lower Baggot street Dublin D02 P593 - Ireland Telephone: +353 1 6971 146 Fax: +353 1 6971 147 Email: ireland@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences AS Address: c/o MerVerdi Munkerudtunet 10 1164 Oslo - Norway Telephone: +47 21 988 882 Email: norge@clinisciences.com Web: https://www.clinisciences.com



Spain Company: CliniSciences Lab Solutions Address: C/ Hermanos del Moral 13

(Bajo E), 28019, Madrid - Spain Telephone: +34 916 750 700 Fax: +34 91 269 40 74 Email: espana@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences Limited Address: 11 Progress Business center, Whittle Parkway, SL1 6DQ Slough- United Kingdom Telephone: +44 (0)1753 866 511 or +44 (0) 330 684 0982 Fax: +44 (0)1753 208 899 Email: uk@clinisciences.com IWeb: https://www.clinisciences.com

Version 07-01-2025



Denmark

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: danmark@clinisciences.com Web: https://www.clinisciences.com



Germany

Company: Biotrend Chemikalien GmbH Address: Wilhelm-Mauser-Str. 41-43, 50827 Köln - Germany Telephone: +49 221 9498 320 Fax: +49 221 9498 325 Email: info@biotrend.com Web: https://www.biotrend.com

Italy

Company: CliniSciences S.r.l Address: Via Maremmana inferiore 378 Roma 00012 Guidonia Montecelio - Italy Telephone: +39 06 94 80 56 71 Fax: +39 06 94 80 00 21 Email: italia@clinisciences.com Web: https://www.clinisciences.com

Poland

Company: CliniSciences sp.Z.o.o. Address: ul. Rotmistrza Witolda Pileckiego 67 lok. 200 - 02-781 Warszawa -Poland Telephone: +48 22 307 0535 Fax: +48 22 307 0532 Email: polska@clinisciences.com Web: https://www.clinisciences.com

Sweden

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: sverige@clinisciences.com Web: https://www.clinisciences.com

USA

Company: Biotrend Chemicals LLC Address: c/o Carr Riggs Ingram, 500 Grand Boulevard, Suite 210 Miramar Beach, FL 32550- USA Telephone: +1 850 650 7790 Fax: +1 850 650 4383 Email: info@biotrend-usa.com Web: https://www.biotrend-usa.com















