Hypherphage

A robust platform to develop recombinant antibodies

Exceptional reproducibility of antibody production

Future-proof supply

Fast production

Controlled conditions

Animal free

Distributed by: CliniSciences Group









The Hyperphage System

An established platform to discover recombinant antibodies, as well as being a powerful tool to improve antibody properties.



PRŒEN



Antibody Library Preparation

The first step is to create antibody variable heavy (VH) and light chain (VL) PCR products. Antibody PCR products are ligated into the pSEX81 surface expression vector (phagemid). This creates a fusion protein containing VH and VL combined with the pIII minor capsid protein derived from the M13 bacteriophage.

Suitable products for this step



- Mouse IgG Library Primer Set, Cat. No. F2010
- Human IgG/IgM Library Primer Set, Cat. No. F2000
- pSEX81 Surface Expression Phagemid Vector, Cat. No. PR3005

Hyperphage Display

Hyperphages carry a deletion in the pIII gene. They are generated by an *E. coli* packaging cell line producing functional pIII. The resulting hyperphages carry functional pIII on their surface but lack the pIII gene in their genome. These hyperphages can be used to infect bacteria with a phagemid library. Due to the resulting display phages carrying several copies of the antibody or peptide on its surface, panning efficiency is increased dramatically.

Suitable products for this step



Quantification

The anti-M13/fd/F1 specifically detects the phage outer surface protein pVIII. Using the Phage Titration ELISA, a standard curve is prepared using a dilution series of phage suspension of known titer. The particle number is estimated in comparision to the standard curve.

Suitable products for this step	
	Phage Titration ELISA, Cat. No. PRPHAGE
	anti-M13/fd/F1 Filamentous Phages mouse monoclonal, B62-FE2, Cat.No. 61097 (available with several conjugates)

Expression of functional recombinant single-chain Fv antibody fragments in *E. coli*

The identified clones of interest can easily be produced by *E. Coli*. This is done by transforming the bacteria with the pOPE101 Expression Vector, which includes the desired antibody fragment. As the expression vector contains a c-Myc and His-tag, the expressed antibody fragment can be purified by metal chelation or affinity purification.

Suitable products for this step



pOPE101 Expression Vector, Cat. No. PR3004

Distributed by: CliniSciences Group



Making gene therapy better. Together.

Our commitment to scientists is to support them in making new scientific discoveries and groundbreaking research within the life sciences and gene therapy. It is our underlying mission, to help progress new therapies safely, quickly and affordably, so together we can provide solutions to people suffering from poor health and affected by disease every day.

PROGEN is made up of a team of experts within AAV and the life sciences and is partnered with gene therapy experts worldwide. We are more than just a manufacturer of antibodies, proteins and ELISA kits. We strive to understand what scientists need, so we can create solutions and ultimately deliver high quality antibodies and exclusive AAV products, which solve your research challenges within academia, biotech and pharma, and ensure PROGEN can continue to be a trusted and reliable partner.

Distributed by: CliniSciences Group

CliniSciences Group

Austria

Company: CliniSciences GmbH Address: Sternwartestrasse 76, A-1180 Wien - Austria Telephone: +43 720 115 580 Fax: +43 720 115 577 Email: oesterreich@clinisciences.com Web: https://www.clinisciences.com

Finland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: suomi@clinisciences.com Web: https://www.clinisciences.com

Iceland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: island@clinisciences.com Web: https://www.clinisciences.com

Netherlands

Company: CliniSciences B.V. Address: Kraijenhoffstraat 137A, 1018RG Amsterdam, - Netherlands Telephone: +31 85 2082 351 Fax: +31 85 2082 353 Email: nederland@clinisciences.com Web: https://www.clinisciences.com

Portugal

Company: Quimigen Unipessoal LDA Address: Rua Almada Negreiros, Lote 5, Loja 14, 2615-275 Alverca Do Ribatejo - Portugal Telephone: +351 30 8808 050 Fax: +351 30 8808 052 Email: info@guimigen.com Web: https://www.quimigen.pt

Switzerland

Company: CliniSciences AG Address: Fracht Ost Flughafen Kloten CH-8058 Zürich - Switzerland Telephone: +41 (044) 805 76 81 Fax: +41 (044) 805 76 75 Email: switzerland@clinisciences.com Web: https://www.clinisciences.com

Belgium

Company: CliniSciences S.R.L Address: Avenue Stalingrad 52, 1000 Brussels - Belgium Telephone: +32 2 31 50 800 Fax: +32 2 31 50 801 Email: belgium@clinisciences.com Web: https://www.clinisciences.com

France

Company: CliniSciences S.A.S Address: 74 Rue des Suisses, 92000 Nanterre- France Telephone: +33 9 77 40 09 09 Fax: +33 9 77 40 10 11 Email: info@clinisciences.com Web: https://www.clinisciences.com

Ireland

Company: CliniSciences Limited Address: Ground Floor, 71 lower Baggot street Dublin D02 P593 - Ireland Telephone: +353 1 6971 146 Fax: +353 1 6971 147 Email: ireland@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences AS Address: c/o MerVerdi Munkerudtunet 10 1164 Oslo - Norway Telephone: +47 21 988 882 Email: norge@clinisciences.com Web: https://www.clinisciences.com

Spain

UK

Company: CliniSciences Lab Solutions Address: C/ Hermanos del Moral 13 (Bajo E), 28019, Madrid - Spain Telephone: +34 916 750 700 Fax: +34 91 269 40 74 Email: espana@clinisciences.com Web: https://www.clinisciences.com



USA

Company: Biotrend Chemicals LLC Address: c/o Carr Riggs Ingram, 500 Grand Boulevard, Suite 210 Miramar Beach, FL 32550- USA Telephone: +1 850 650 7790 Fax: +1 850 650 4383 Email: info@biotrend-usa.com Web: https://www.biotrend-usa.com



Denmark

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: danmark@clinisciences.com Web: https://www.clinisciences.com



Germany

Company: Biotrend Chemikalien GmbH Address: Wilhelm-Mauser-Str. 41-43, 50827 Köln - Germany Telephone: +49 221 9498 320 Fax: +49 221 9498 325 Email: info@biotrend.com Web: https://www.biotrend.com



Italy

Company: CliniSciences S.r.I Address: Via Maremmana inferiore 378 Roma 00012 Guidonia Montecelio - Italy Telephone: +39 06 94 80 56 71 Fax: +39 06 94 80 00 21 Email: italia@clinisciences.com Web: https://www.clinisciences.com

Poland

Company: CliniSciences sp.Z.o.o. Address: ul. Rotmistrza Witolda Pileckiego 67 lok. 200 - 02-781 Warszawa -Poland Telephone: +48 22 307 0535 Fax: +48 22 307 0532 Email: polska@clinisciences.com Web: https://www.clinisciences.com

Sweden

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064 Email: sverige@clinisciences.com

Web: https://www.clinisciences.com









Version 07-01-2025