

# **IMMUNOHISTOCHEMISTRY BROCHURE**

Distributed by:



Distributed by:



## Introduction

Immunohistochemistry (IHC) is a method for detecting specific antigens present on tissues through the use of purified antibodies and chromogen detection systems. The IHC staining process can seem complex and requires many different standardized reagents and ancillaries. As experts in IHC, ScyTek Laboratories can provide the optimal set of reagents for your specific needs. We offer traditional methods, such as streptavidin-biotin systems, as well as more sensitive and modern options, such as secondary polymers.

Distributed by:



ScyTek Laboratories offers over **500** antibodies for immunohistochemistry. We're confident we can provide you with the right antibody for your experiment. Offerings include:

- Mouse or Rabbit. Polyclonal and Monoclonal.
- Both pre-dilute/ready-to-use and concentrated format
- Concentrated sizes available as 0.1ml or 1ml
- Ready-to-use sizes available up to 1L
- Validation data, technical support, and guarantee of performance




Availability Sizes		Format	
"A" Antibodies RTU: 2mls, 7mls, 25mls Concentrate: 0.1ml, 1ml		RTU and Concentrate	
"RA" Antibodies (Item#: RA00XXX)	0.5 mls, 1 ml	Concentrate only	


RTU: Ready-to-use, X: Any number


# OUR MOST PROMINENT PRIMARY ANTIBODIES

<b>Antibody Description</b>	Item #
CD20, B-Cell; Clone L26	A00003
CD3, T-Cell; Polyclonal (Rabbit)	A00030
Cytokeratin 7; Clone OV-TL12/30	A00142
S-100; Polyclonal (Rabbit)	A00036
Epithelial Membrane Antigen; Clone E29	A00008
Ki-67 Antigen; Polyclonal	A00095
CD45, Leucocyte Common Antigen(LCA); Clones PD7/26 & 2B11	A00017
Desmin; Clone D33	A00007
Actin, Alpha-Smooth Muscle; Clone1A4	A00002
c-erbB-2 Oncoprotein; Clone SP3	A00100
Cytokeratin, Pan; Clones AE1 & AE3	A00152
MIC2 Gene Protein, CD99; Clone HO36-1.1	A00044



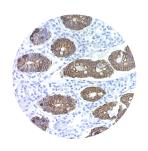


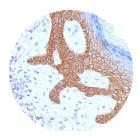

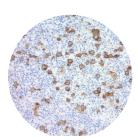


Research Use Only Antibodies	Item #
CD56 / NCAM1 / NKH1 (Neuronal Cell Marker); Clone 123A8 (56C04)	RA0238
p53 Tumor Suppressor Protein; Clone BP53-12 & DO-7	RA0321
MUC2 (Mucin 2); Clone CCP58	RA0219
Caldesmon, HMW (h-Caldesmon) (Smooth Muscle Marker); Clone h-CALD	RA0028
CDX2 (GI Epithelial Marker); Clone CDX2/1690	RA0480
Chromogranin A / CHGA (Neuroendocrine Marker); Clone LK2H10, PHE5 & CGA/414	RA0095
PDCD1 / PD1 / CD279 (Programmed Cell Death 1); Clone NAT105	RA0254
CD117/c-Kit (Marker for Gastrointestinal Stromal Tumors); Clone C117/370	RA0168
Ksp-Cadherin (Kidney-Specific Cadherin) / CDH16; Clone CDH16/1532R	RA0470
Complement 4d (C4d) (Acute Humoral Rejection Marker); CloneC4D204	RA0025

Distributed by:

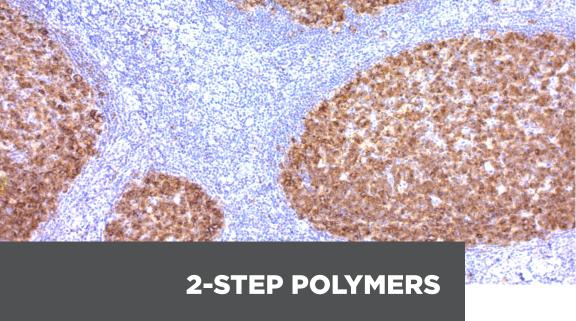

Save time and energy spent on 3-step IHC processes by using modern 2-step Polymer reagents. Our polymer detection reagents are more sensitive than traditional methods and come in both rabbit and mouse specific formats. Offerings include:

- Mouse, Rabbit, and Polyvalent
- Both HRP and Alk-Phos conjugation
- Full IHC Kits containing Polymers
- Wide range of volumes and kit configurations offered

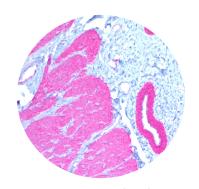



Distributed by:
CliniSciences Group

## Single Polymer Solutions

Description	Item Code:
CRF Anti-Polyvalent HRP Polymer	ABZ
PolyTek Anti-Mouse Polymerized Alk-Phos	PAT
PolyTek, Anti-Mouse Polymerized HRP	PAM
PolyTek, Anti-Rabbit Polymerized HRP	PAR
PolyTek, Anti-Rabbit Polymerized Alk-Phos	PAL

Kits Offered:	Item #:	# of Slides
CRF Anti-Polyvalent HRP Polymer (DAB)	CPH080	80
Stain Kit		
Contents: Peroxide Block 8ml, Super Block 8ml,		
HRP Polymer 8ml, DAB Substrate 7 x 5ml, DAB		
Chromogen 3ml.		
CRF Anti-Polyvalent HRP Polymer (DAB) Lab	CPP125	1,250
Pack	CPP500	5,000
Contents: Super Block, Anti-Polyvalent HRP Poly-	CPP999	10,000
mer, DAB Substrate, DAB Chromogen		
D. I. T. I. VIDD. I I	PIDOGO	
PolyTek HRP Anti-Rabbit Polymerized Imaging	PIP080	70
System		
Contents: Peroxide Block for Image Analysis - 1 x		
8ml - ADA008		
Super Block - 1 x 8ml - AAA008 CRF Anti-Rabbit HRP - 1 x 8ml PAR008		
DAB Chromogen - $1 \times 3ml$ - $ACB003$		
DAB Substrate (High Contrast) - 8 x 5ml - ACU005		
DAD Substrate (High Contrast) - 6 x 5mt - ACC005		
PolyTek Anti-Mouse (DAB) Polymerized HRP	PIR080	70
Imaging System	1111000	, 0
Contents: Peroxide Block for Image Analysis - 1 x		
8ml - ADA008		
Super Block - 1 x 8ml - AAA008		
CRF Anti-Mouse HRP - 1 x 8ml - PAM008		
DAB Chromogen - 1 x 3ml - ACB003		
DAB Substrate (High Contrast) - 8 x 5ml - ACU005		
Distributed by:		

# STREPTAVIDIN/BIOTIN SYSTEMS

These reagents are for those who prefer the more traditional and economical method of staining. There are several choices of species and concentrations of biotinylated secondaries for your IHC procedure. Offerings include:

 Anti-Mouse, Anti-Rabbit, Anti-Human, and Anti-Goat for secondary antibodies

- Combined specificity reagents known as 'Polyvalent'
- Alk-Phos and HRP enzymes

Differing formats of these systems are offered at respective prices, allowing users to choose more economical options while still achieving excellent results through slightly longer staining times.

	(standard	(higher	(highest
	sensitivity)	sensitivity)	sensitivity)
Suggested Staining Time (Minutes):	30	20	10

# Steptavidin/Biotin Kit Examples:

	Description	Item #	Slides
	UltraTek HRP Anti-Polyvalent (DAB)	AMF080	70
	Staining System		
	UltraTek HRP Anti-Polyvalent Staining	AFN600	500
	System		
	UltraTek HRP Anti-Polyvalent Lab Pack	UHP125	1,250
	UltraTek Complete HRP Anti-Polyvalent	UCS125	1,250
	(DAB) Staining System	. =====	
	UltraTek Alk-Phos Anti-Mouse (Permanent Red) Staining System	AET080	70
	UltraTek Alk-Phos Anti-Mouse Lab Pack	UAM125	1,250
	UltraTek Alk-Phos Anti-Rabbit (Permanent	AME080	70
	Red) Staining System		
	UltraTek Alk-Phos Anti-Rabbit Staining	AFM600	500
	System		
	UltraTek HRP Anti-Polyvalent (AEC)	AMG080	70
	Staining System		
	SensiTek HRP Anti-Polyvalent Lab Pack	SHP125	1,250
	SensiTek HRP Anti-Polyvalent (DAB)	AEM080	70
	Staining System		
	SensiTek Alk-Phos Anti-Polyvalent (Per-	AES080	70
"园	manent Red) Staining System		
10	SensiTek Alk-Phos Anti-Rabbit Staining	AFF600	500
	System		
	SensiTek HRP Anti-Mouse Staining Sys-	AFA600	500
	tem		
	SensiTek HRP Anti-Rabbit (AEC) Staining	AEJ080	70
	System		
	EconoTek HRP Anti-Polyvalent (DAB)	AEX080	70
	EconoTek Alk-Phos Anti-Polyvalent Lab	EAP125	1,250
	Pack		1.5-
	EconoTek HRP Anti-Polyvalent Lab Pack	EHP125	1,250

A substantial amount of individual reagents and kit sizes/combinations are available. Please visit website or contact us for entire list of offerings.



One key strength is our ability to provide most IHC reagents in several different combined kit formats at several different sizes. This flexibility, along with offering nearly any size of individual reagent, allows us to cater to your specific needs and requirements. Thousands of liters per year of ScyTek's premium DAB chromogen/substrate solutions are used by clinicians and researchers around the world. Offerings include:

- Kits containing most reagents needed for IHC link, label, chromogen/substrate, and counterstain
- Bulk chromogen/substrate Kits
- Standard DAB, Permanent Red, and AEC options
- BCIP NBT and other colors
- Standard and high sensitivity options



Distributed by:

# Popular DAB Kits containing substrate and chromogen

	Description	Item #	# of Slides*
lard	DAB Chromogen/Substrate	ACH500	500
Standa	DAB Chromogen/Substrate Bulk pack	ACK500	5,000
St		ACK999	1,000
	DAB Chromogen/Substrate Bulk Pack	ACV500	5,000
ligh	(High Contrast)	ACV999	
iI	DAB Chromogen/Substrate Kit (High	ACT500	500
	Contrast)		

<sup>\*</sup>Slides stained based on a 100µl basis per slide

Description	Item #:	# of Slides
Permanent Red Kit (For Alk-Phos)	PRD-15	150
	PRD500	500
AEC Chromogen/Substrate Kit	ACJ500	5,000
	ACJ999	10,000


When ordering Chromogen and Detection system kits in bulk sizes ('Bulk Kit', 'Lab Packs', 'Bulk Pack') a corresponding kit must be ordered to receive proportionate volume of reagents. Below are some examples:

If you order:	You may also need:
AFN600	ACH500/ACT500
UHP	ACK/ACV

We also offer an IHC kit that contains the hematoxylin counterstain:

UltraTek Complete System; Anti-Polyvalent, HRP, DAB	Item #:
Contents: Peroxide block. Super block, Link, Label,	UCS015
Chromogen, Hematoxylin, Bluing Reagent	UCS125




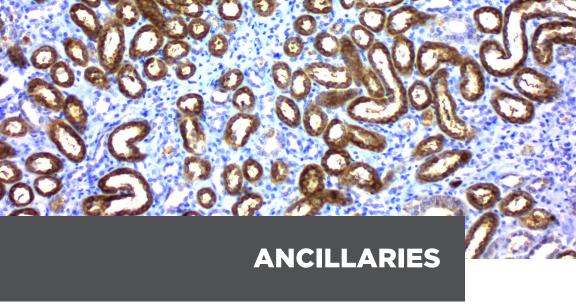

IHC is a complex process that requires many other dependable reagents for a successful stain. We offer nearly all of those reagents.

- Antigen retrieval EDTA, tris, citrate, high ph, low ph
- Enzymatic retrievals trypsin, pepsin
- Wash Buffers tris, phosphate, economical concentrated solutions (10X to 25X), powdered
- Antibody diluents, stabilizing diluents
- Blocking solutions peroxidase, protein, mouse, rabbit, and human
- Mounting media
- Formalin fixatives
- Marking pens, PAP pens
- Counterstains



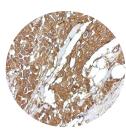

# Popular Ancillaries:

<b>Blocking Reagents</b>	Item code:
Peroxide Block	ACA
Super Block	AAA
Biotin Blocking Kit for Image Analysis	BBK
Human-to-Human Blocking Reagent	НТН
Rabbit-to-Rabbit Blocking Reagent	RTR

# Popular Ancillaries (cont'd):


IHC Wash Buffers	Item code:
Phosphate Buffered Saline plus Tween 20 (20x) pH 7.6	PBT
Tris Buffered Saline plus Tween 20 (20x Concentrate) pH 7.4	TBT
Phosphate Buffered Saline plus Tween 20 (10x) pH 7.4	PBE
Tris Buffered Saline plus Tween 20 (10x) pH 7.5	TBE
Wash buffers without Tween are also offered.	

Retrieval Solutions	Item code:
Pepsin, Stabilized Solution	PSS
Trypsin Enzymatic Digestion Kit	TSS
Tris-EDTA HIER Solution (10x) pH 9.0	TES
Citrate Buffer (10x) pH 6.0	CBB
EDTA - Saline Buffer (10X Concentrate); pH 8.0	ETA
Citrate Plus (10x) HIER Solution	CPL

Primary Antibody Diluent	Item code:
Primary Antibody Diluent (Phosphate, Green)	APG
Primary Antibody Diluent (Tris, Green)	ATG
Colorless diluent is offered as well.	

Counterstaining	Item code:
Hematoxylin, Mayer's (Lillie's Modification)	HMM
Bluing Reagent	BRT
Hematoxylin and Eosin Stain Kit	HAE
Nuclear Fast Red Solution	NFS

Other	Item code:
DAB Away Kit	ACP001
DAB (Post) Enhancing Solution	DES

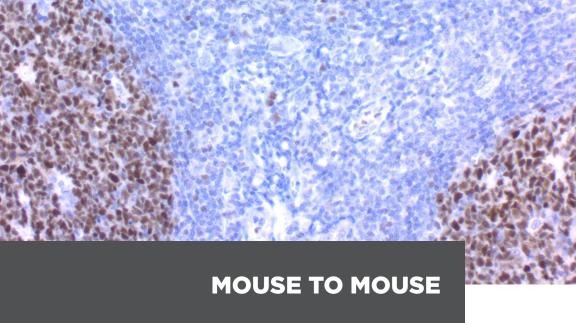

ScyTek offers optimized blocking reagents of endogenous Ig, allowing for staining on tissues that are the same species as the primary antibody. Our mouse to mouse kit has been a very popular and effective option for IHC staining on mouse tissue. Offerings include:

- Entire kits containing detection systems and chromogen for either HRP or Alk phos
- Mouse, Rabbit, and Human Blocking reagents



Distributed by:
CliniSciences Group

# Example of Kits Offered:

Item Code

Mouse to Mouse HRP (DAB) Staining System

MTM001

Contents:

Peroxide Block - 1 x 8ml

Super Block - 1 x 8ml

Mouse Block - 1 x 8ml

Anti-Polyvalent - 1 x 8ml

HRP - 1 x 8ml

DAB Concentrate - 1 x 3ml

DAB Substrate Buffer - 6 x 5ml

MTM004

Mouse to Mouse Alk-Phos (Permanent

Red) Staining System

Contents:

Super Block - 1 x 8ml

Mouse Block- 1 x 8ml

Anti-Polyvalent - 1 x 8ml

Alk-Phos - 1 x 8ml

Permanent Red - 1 x 2ml

Permanent Red Buffer - 7 x 5ml

Mouse to Mouse HRP (AEC) Staining System

MTM002

Contents:

Peroxide Block - 1 x 8ml

Super Block - 1 x 8ml

Mouse Block - 1 x 8ml

Anti-Polyvalent - 1 x 8 ml

HRP - 1 x 8ml

AEC Chromogen - 1 x 3ml

AEC Substrate Buffer - 6 x 5ml

Distributed by:

All of ScyTek's reagents for IHC are made with high quality proteins and purification methods. All other raw materials are of ACS grade or higher and water used is verified to conform to industry standards. All batches of reagents and kits undergo a rigorous quality control procedure before release. Strict validation procedures of each reagent and kit have been developed by our scientists to ensure that results are in line with published literature and/or expected results stated on the Instructions For Use (IFU).



## Technical Support:

If any results appear inconsistent with what should be expected, please contact by phone or email ÿour local distributor

Distributed by:

# CliniSciences Group

### Austria

Company: CliniSciences GmbH Address: Sternwartestrasse 76, A-1180

Wien - Austria

Telephone: +43 720 115 580 Fax: +43 720 115 577

Email: oesterreich@clinisciences.com Web: https://www.clinisciences.com



### Belgium

Company: CliniSciences S.R.L Address: Avenue Stalingrad 52, 1000

Brussels - Belgium Telephone: +32 2 31 50 800 Fax: +32 2 31 50 801

Email: belgium@clinisciences.com Web: https://www.clinisciences.com



### Denmark

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: danmark@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: suomi@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences S.A.S. Address: 74 Rue des Suisses, 92000

Nanterre- France

Telephone: +33 9 77 40 09 09 Fax: +33 9 77 40 10 11 Email: info@clinisciences.com

Web: https://www.clinisciences.com



Company: Biotrend Chemikalien GmbH Address: Wilhelm-Mauser-Str. 41-43. 50827 Köln - Germany Telephone: +49 221 9498 320

Fax: +49 221 9498 325 Email: info@biotrend.com Web: https://www.biotrend.com

### Iceland

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: island@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences Limited Address: Ground Floor, 71 lower Baggot street Dublin D02 P593 - Ireland

Telephone: +353 1 6971 146 Fax: +353 1 6971 147

Email: ireland@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences S.r.l Address: Via Maremmana inferiore 378 Roma 00012 Guidonia Montecelio - Italy Telephone: +39 06 94 80 56 71 Fax: +39 06 94 80 00 21

Email: italia@clinisciences.com Web: https://www.clinisciences.com

### Netherlands

Company: CliniSciences B.V. Address: Kraijenhoffstraat 137A 1018RG Amsterdam, Netherlands Telephone: +31 85 2082 351 Fax: +31 85 2082 353

Email: nederland@clinisciences.com Web: https://www.clinisciences.com



### Norway

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: norge@clinisciences.com Web: https://www.clinisciences.com



Company: CliniSciences sp.Z.o.o. Address: ul. Rotmistrza Witolda Pileckiego 67 lok. 200 - 02-781 Warszawa -Poland Telephone: +48 22 307 0535

Fax: +48 22 307 0532 Email: polska@clinisciences.com

Web: https://www.clinisciences.com



Company: Quimigen Unipessoal LDA Address: Rua Almada Negreiros, Lote 5, Loia 14, 2615-275 Alverca Do Ribatejo - Portugal Telephone: +351 30 8808 050

Fax: +351 30 8808 052 Email: info@quimigen.com Web: https://www.quimigen.pt



### Spain

Company: CliniSciences Lab Solutions Address: C/ Hermanos del Moral 13 (Bajo E), 28019, Madrid - Spain Telephone: +34 91 269 40 65 Fax: +34 91 269 40 74

Email: espana@clinisciences.com Web: https://www.clinisciences.com

Company: CliniSciences Limited



### Sweden

Company: CliniSciences ApS Address: Oesterbrogade 226, st. 1, Copenhagen, 2100 - Denmark Telephone: +45 89 888 349 Fax: +45 89 884 064

Email: sverige@clinisciences.com Web: https://www.clinisciences.com



### Switzerland

Company: CliniSciences Limited Address: Marktgasse 18 8302 Kloten -Switzerland

Telephone: +41 (044) 805 76 81 Fax: +41 (044) 805 76 75

Email: switzerland@clinisciences.com Web: https://www.clinisciences.com



Address: 11 Progress Business center, Whittle Parkway, SL1 6DQ Slough- United Kingdom Telephone: +44 (0)1753 866 511

or +44 (0) 330 684 0982 Fax: +44 (0)1753 208 899 Email: uk@clinisciences.com Web: https://www.clinisciences.com



Company: Biotrend Chemicals LLC Address: c/o Carr Riggs Ingram, 500 Grand Boulevard, Suite 210 Miramar Beach, FL 32550- USA Telephone: +1 850 650 7790 Fax: +1 850 650 4383 Email: info@biotrend-usa.com

Web: https://www.biotrend-usa.com

